

Accessing Frozen Room Leica LV1 via GotoAssist:

- **Local user (at SHY)** logs in as LCA-LV1-SHY\aperio and Sc@nscope1234
- **Local user (at SHY)** launches the LV1 Console (Scanner Application)
-
- **Local user (at SHY)** loads the Slides
- **Remote user** visits <http://app.gotoassist.com>
- **Remote user** will use UPMC email address as the Username and the password they have created at GotoAssist (*Not the UPMC Password*)
- **Remote user** will be on the Dashboard
- **Remote user** will Click the *BLUE Connect* button next to LCA-LV1-SHY PC name
- **Remote user** logs in to the same desktop as LCA-LV1-SHY\aperio and Sc@nscope1234
- **Remote user** will be on the same Desktop profile as the Local user
- **Remote user** will end the GotoAssist session by selecting End Session (Top left corner, pull-down menu and select End Session)

